

Dip Sampler, BÜRKLE

Dip sampler for liquids

Easy sampling of liquid samples. Good view through transparent tubes. Upper end with stopcock, lower end with ball valve.

- PP, transparent
- Dip sampler 100455200 is continuously extractable from 100 cm to 200 cm
- All-layer sample and target-point sample possible

Function

All-layer sample 1. Slowly insert dip sampler in liquid with opened stopcock .
2. Close valve. 3. Remove dip sampler and take sample.

Target-point sample 1. Insert dip sampler in target point with closed stopcock.
2. Open stopcock - liquid pours in. 3. Close stopcock. 4. Remove dip sampler, take sample.

	Volume ml	Immersion depth cm	Ø mm	Order No.
①	140	50	25	10 04 55050
②	280	100	25	10 04 55100
③	500	100 – 200	25	10 04 55200



WWR4Q050

